FFT and Spectrum Analyzer

Objective:

- Understand the frequency domain and some of it's uses
- Understand the Discrete Fourier Transform

Preparation: (i) PC and data acquisition system (ii)FFT software

Background:

A signal can be observed from two different domains:

- 1. The time domain
- 2. The frequency domain

A spectrometer, e.g. a prism, splits light into its component colors which provides its spectral content. This is similar to the trace on a spectrum analyzer, where the abscissa is the frequency and the ordinate is the signals amplitude at that frequency. We have been familiar with the *time domain* signal such as the trace on an oscilloscope where the ordinate is the signals amplitude, and the abscissa is the time variable.

Any signal can be fully described in either of these domains and they are related by a tool called the *Fourier transform*. The Discrete Fourier Transform (DFT) is used to produce frequency analysis of discrete non-periodic signals.

It is usually employed to transform data which may arise from either an actual continuous time process or perhaps a discrete time process which is being analysed from a continuous time system approach. The DFT is usually used to approximate the Fourier transform of a continuous time process, and it is necessary to understand some of the limitations inherent in this approach.

There are three possible phenomena that result in errors between the computed and the desired transform. These three phenomena are (a) *aliasing*, (b) *leakage*, and (c) *the picket-fence effect*.

(a) *Aliasing*. The only solution to the aliasing problem is to ensure that the sampling rate is high enough to avoid any spectral overlap, or to use an *anti-aliasing filter*.

(b) *Leakage*. This problem arises because of the practical requirement of the signal to be a finite interval. The process of terminating the signal after a finite number of terms is equivalent to multiplying the signal by a *window function*. The net effect is a

distortion of the spectrum. There is a spreading or leakage of the spectral components away from the correct frequency, resulting in an undesirable modification of the total spectrum.


(c) *Picket-Fence Effect.* This effect is produced by the inability of the DFT to observe the spectrum as a continuous function, since computation of the spectrum is limited to integer multiples of the fundamental frequency F (reciprocal of the sample length). Observation of the spectrum with the DFT is analogous to looking at it through a sort of "picket-fence," since we can observe the exact behavior only at discrete points. The major peak of a particular component could lie between two of the discrete transform lines, and the peak of this component might not be detected without some addition processing.

A fast Fourier transform (FFT) is an efficient algorithm to compute the DFT and its inverse. It reduces the number of computations needed for N points from $2N^2$ to $2N\log_2 N$.


Procedure:


- 1. Turn on the PC. Run the program "Sigview."
- 2. Follow the procedure given below to be familiar with the software.
- 3. Observe the results carefully and pay attention to the *sampling* (sampling rate) and *truncation* (data length) effect.
- 4. Generate your own signals and perform FFT to observe their spectra such that you can complete your report.
- 5. Use sinus and step functions only.


🕵 Sigview		- 8 🛛
<u>File</u> <u>Data acquisition</u> <u>Edit</u> <u>Play & navigate</u>	Signal tools <u>3D</u> tools <u>System control</u> <u>W</u> indow <u>H</u> elp	
🖻 🙀 🖬 🗃 🎆 🛼 🛼 🛛 K	Spectral analysis defaults	止業■☆ 鮎上手雪♪ 図醸 !
	Iff. EFT spectrum analysis Ctrl+F Iff. Filter Smoothing Remove linear trend Remove peaks Peak hold Averager	
	Ime FFT Ctil+T Spectrogram Track changes as 3D graphics	
	Use custom tool	
	Instruments and markers Log instrument values	
	Signal generator	
	📰 Signal <u>c</u> alculator	
Generate standard wave signals		
🛃 開始 🥂 🧟 Sigview		🖮 🕥 🛃 🍐 上午 09:26

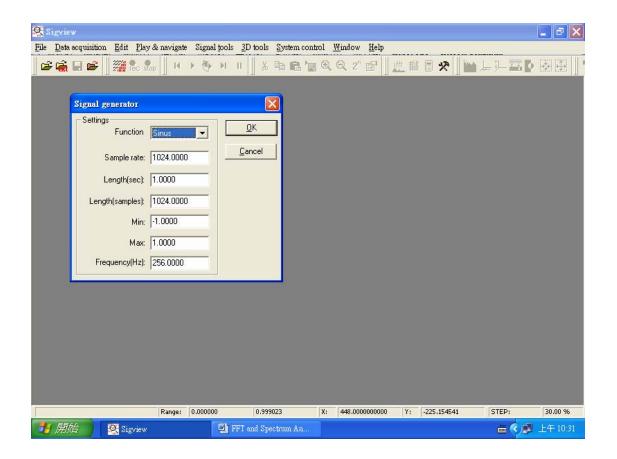



🕵 Sigview			- 7 🗙
File Data acquisition Edit Play & navigate	Signal tools 3D tools System control		
🛛 🖙 🚒 🖬 📽 🖉 🎆 🗞 🗞 📗 M	Spectral analysis defaults	🖆 📗 🏨 🖩 🛠 📗 🖕 斗 🧰 🖡	
421CEN (Sin) 08 08 04 04 02 04 02 04 02 04 02 04 02 04 02 04 05 04 05 04 05 05 06 01 02 02 03 04 05 05 01 02 02 03 04 05 05 02 03 04 05 04 05 04 05 04 05 04 05 04 05 04 05	Filter Smoothing Remove linear trend Remove peaks Peak hold Averager Time FFT Spectrogram Track changes as 3D graphics Use custom tool Instruments and markers Log instrument values Signal generator Signal calculator Signal calculator	Cttl+T	
Spectrum of visible part of the signal			
🛃 開始 🧟 Sigview	🖳 FFT and Spectrum An		上午 10:32

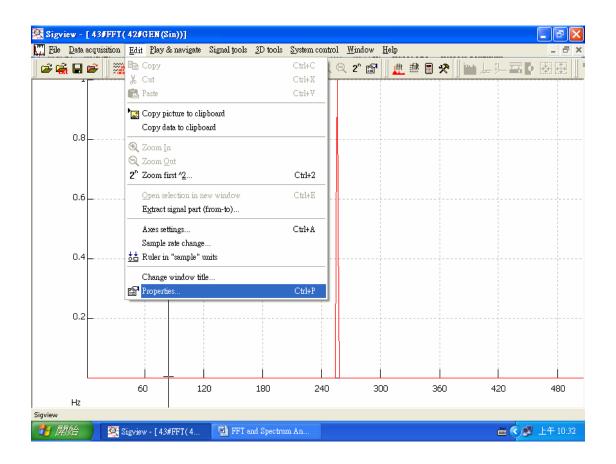


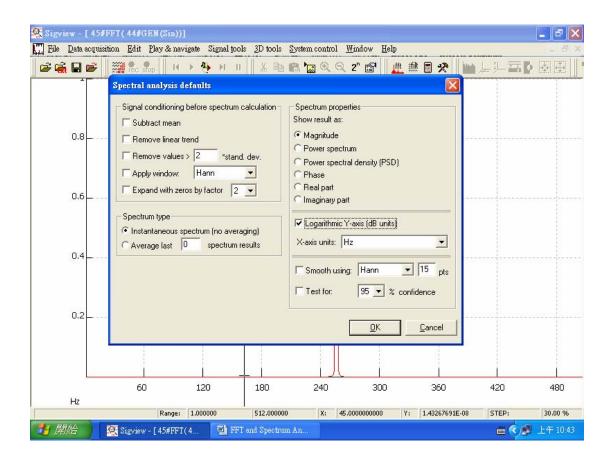


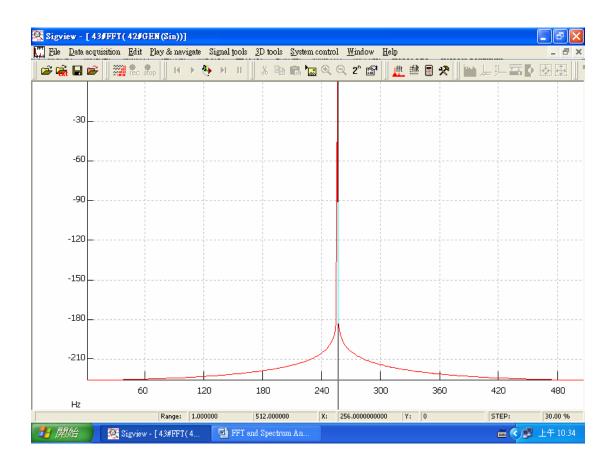


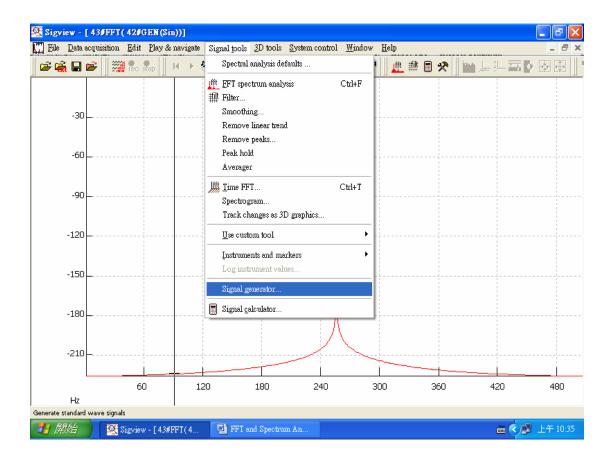


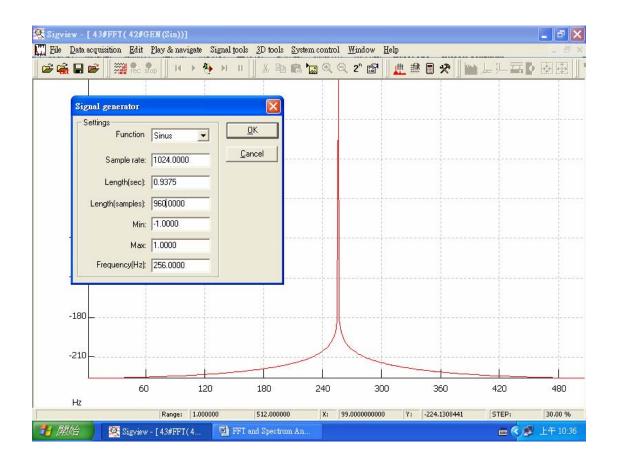


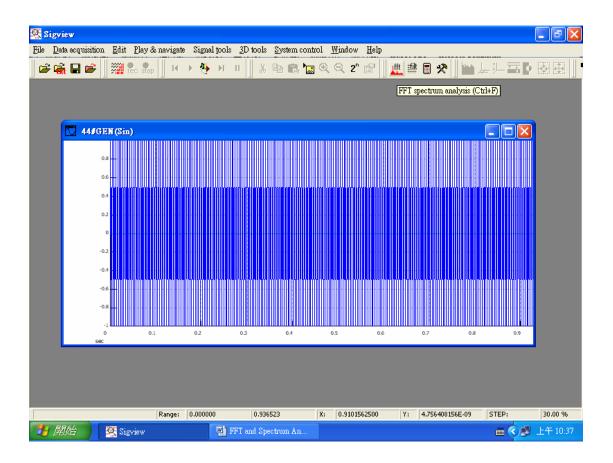


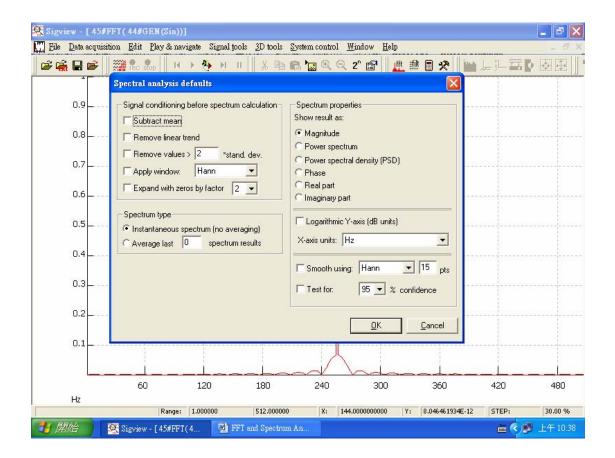


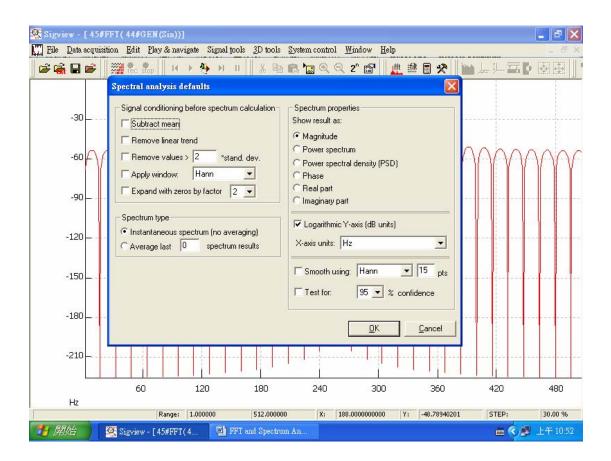


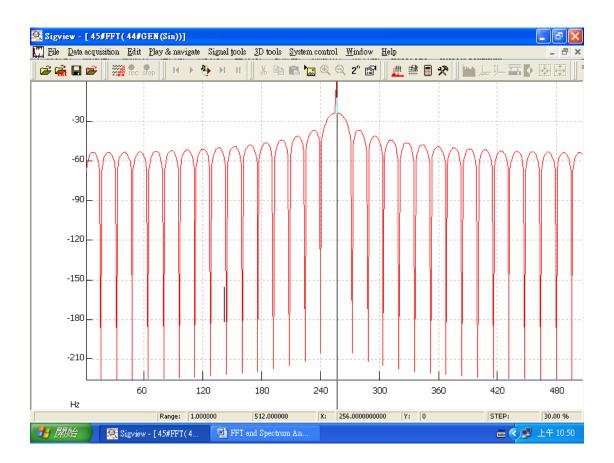


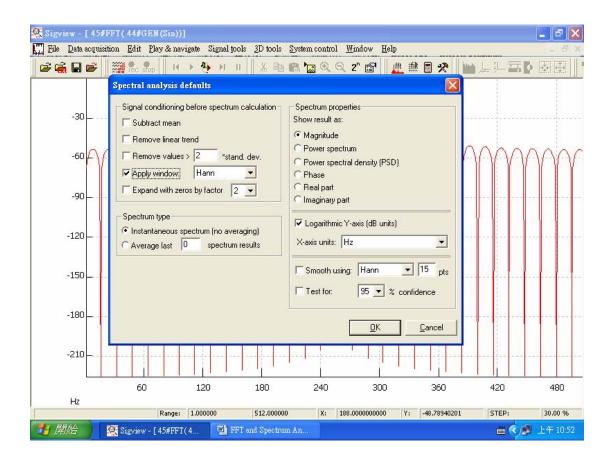


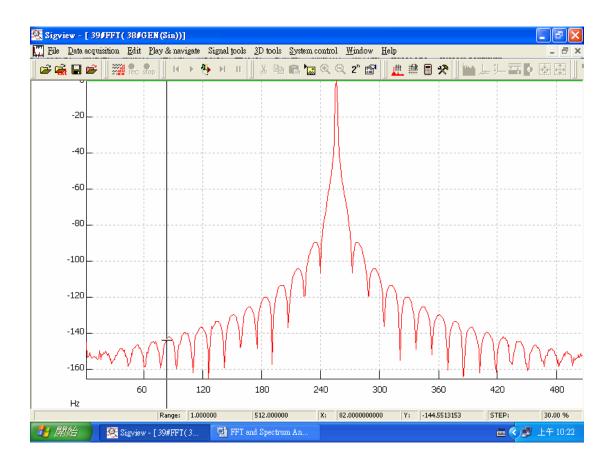


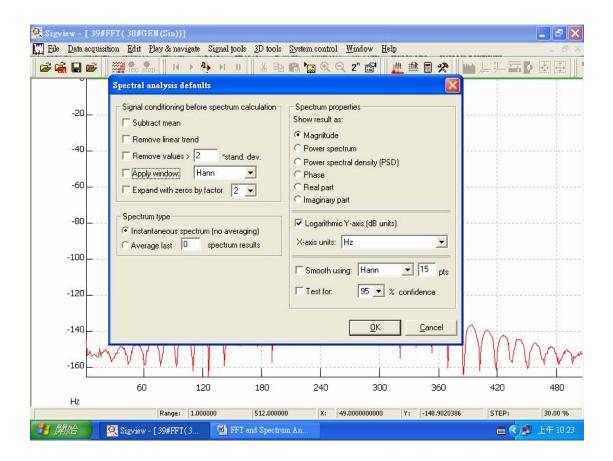


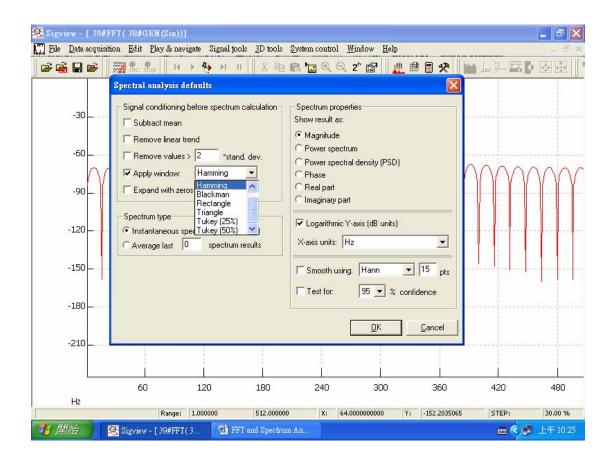


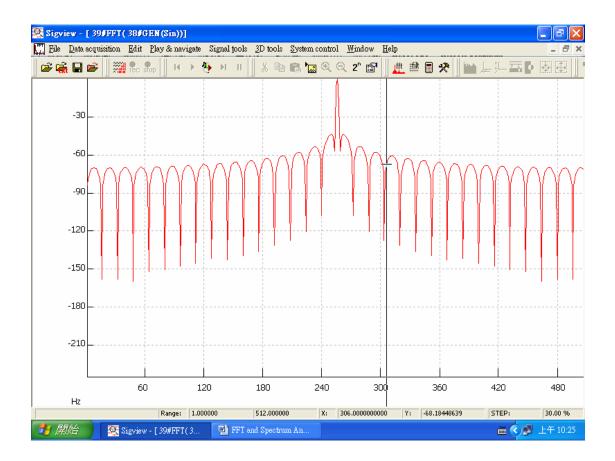


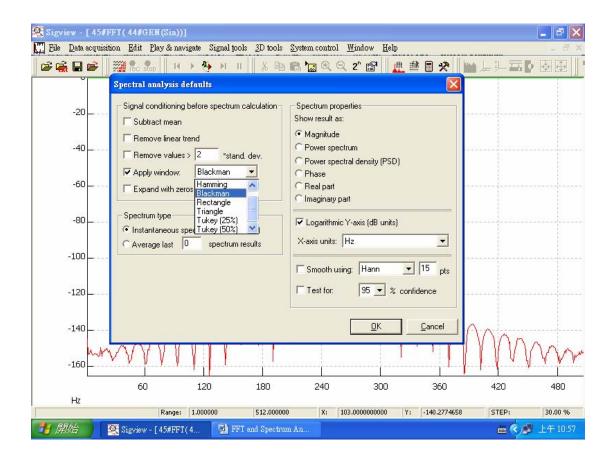


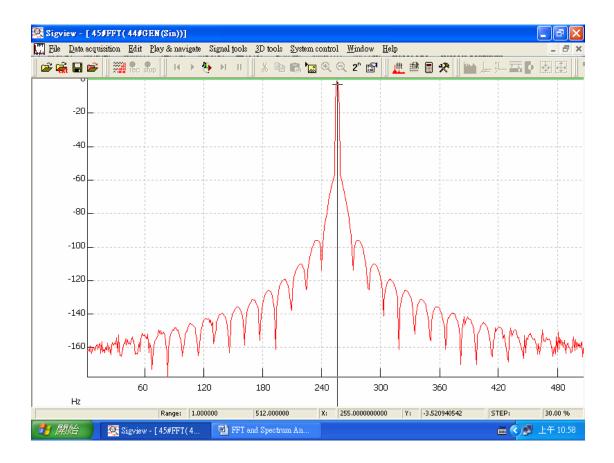


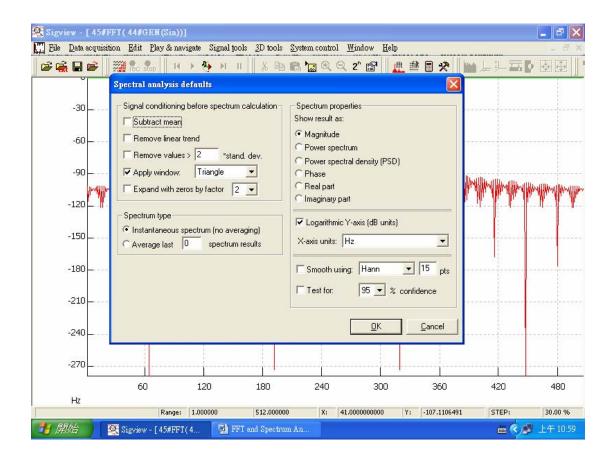


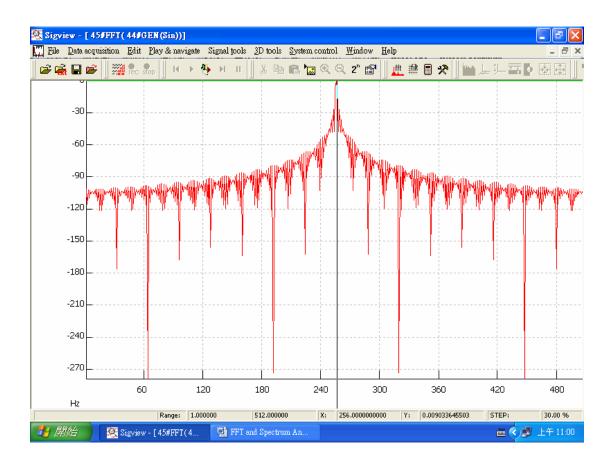


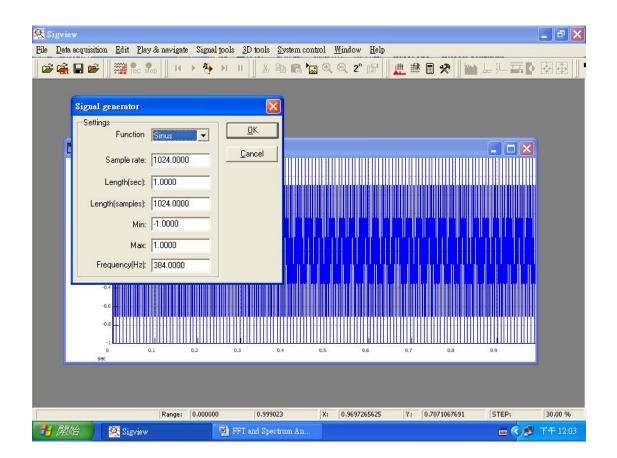


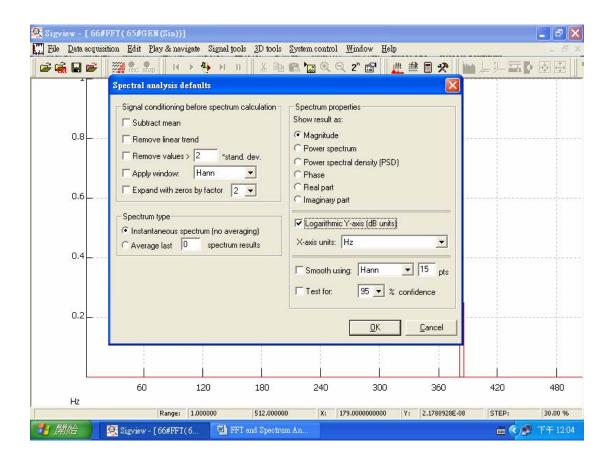


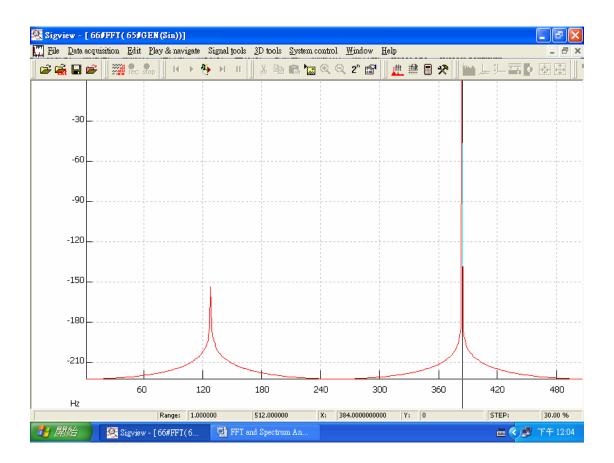


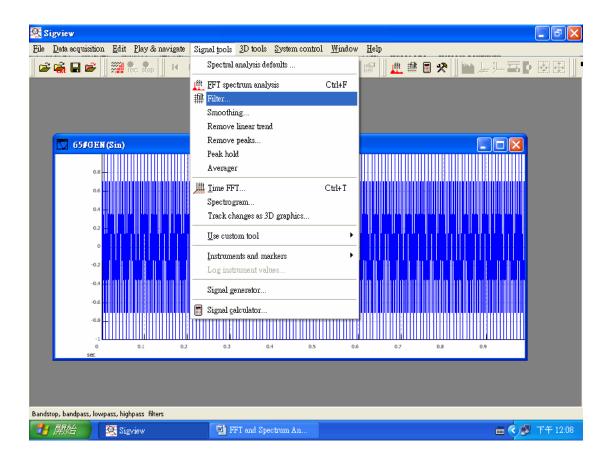


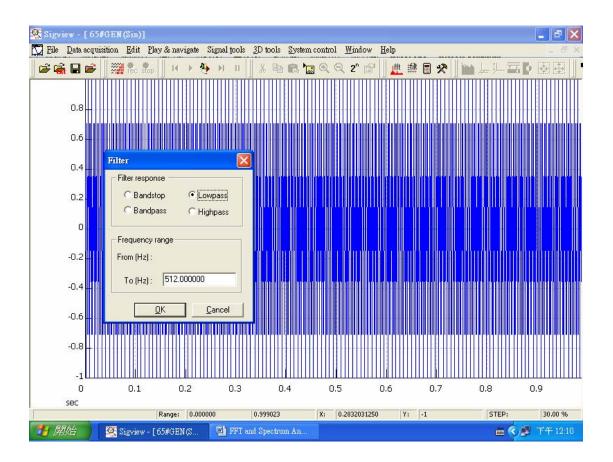

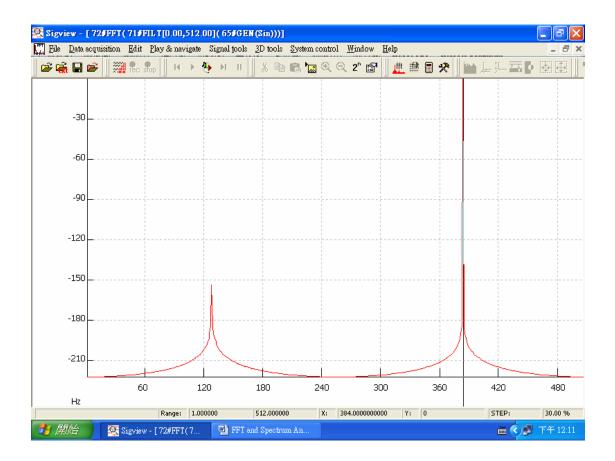

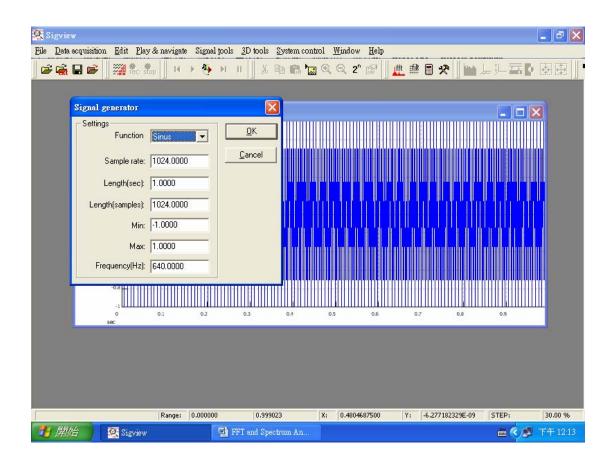


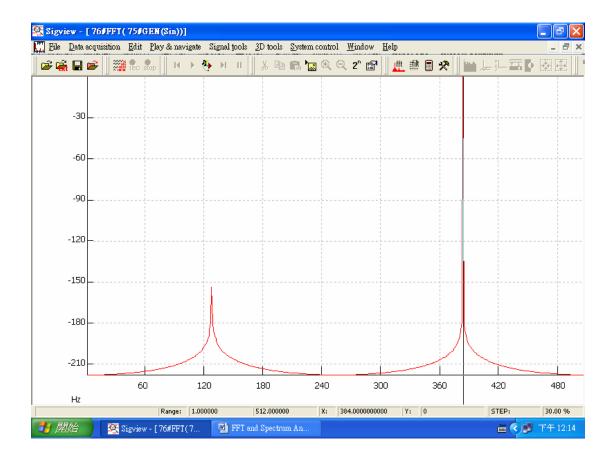


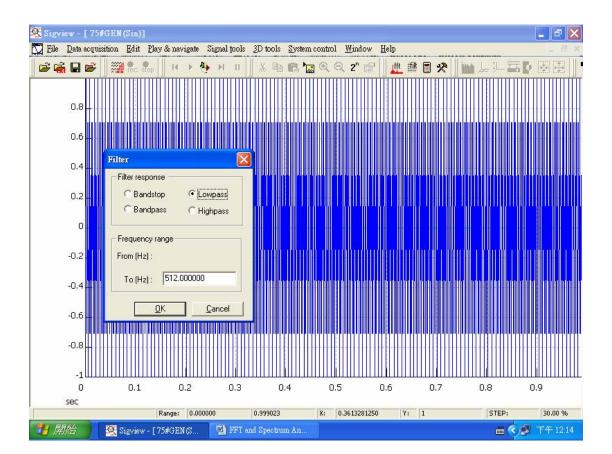


🕵 Sigview					∎₽⊠
<u>File</u> <u>D</u> ata acquisition <u>E</u> dit <u>P</u> lay & navigate	Signal tools <u>3D tools</u> System contro	l <u>W</u> indow <u>H</u> elp)		
🖙 🙀 🖬 🗃 🦷 🕅 🗰 🗛	Spectral analysis defaults	r 🖉	<u>#</u> #E) 🛠 🛛 🔛 🔎	二百 図題 !!!
	Iff. EFT spectrum analysis ## Filter Smoothing Remove linear trend Remove peaks Peak hold Averager	Cttl+F	-	12	
	Jime FFT Spectrogram Track changes as 3D graphics	Ctrl+T			
	Use custom tool	•			
	Instruments and markers Log instrument values	•			
	Signal generator				
	Signal <u>c</u> alculator				
Tapering, smoothing,					
🛂 開始 🔤 🖬 FFT and Spectrum Ar	🔯 Sigview				🛗 🔇 🦻 下午 04:28

Sigview File Data acquisition	Elia Dine Summer Completela 2014-1- Co	star sector Widday The	
File Data acquisition	Edit Play & navigate Signal tools 3D tools Sy		
	Spectral analysis defaults		
	Signal conditioning before spectrum calculation Subtract mean Remove linear trend Remove values > 2 *stand. dev. Apply window: Hann • Expand with zeros by factor 2 • Spectrum type Instantaneous spectrum (no averaging) Average last 0 spectrum results	Spectrum properties Show result as: Magnitude Power spectrum Power spectral density (PSD) Phase Real part Magnithmic Y-axis (dB units) X-axis units: Hz Smooth using: Hann 15 pts Test for: 95 2 % confidence	
		<u> </u>	
	Range: 0.000000 0.999023	X: 0.9638671875 Y: 5.036975281E-09	STEP: 30.00 %
🛃 開始 🔰	🐏 FFT and Spectrum An 🦉 Sigview		🖮 📢 🦻 下午 04:29

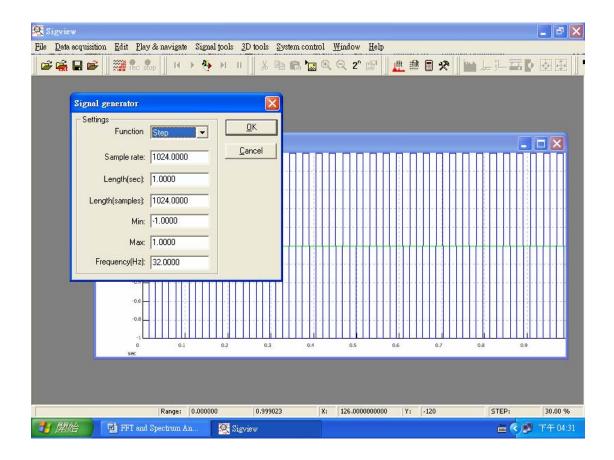

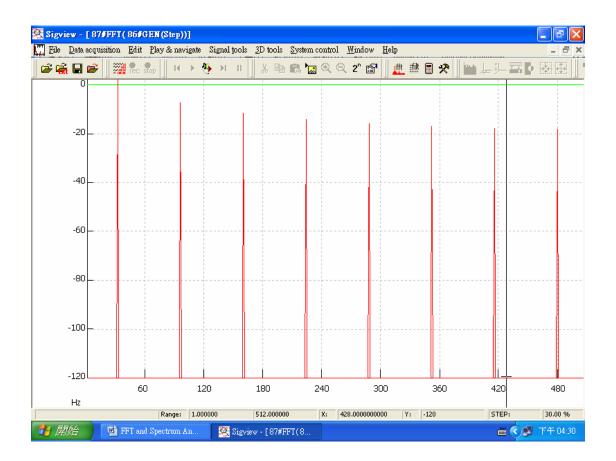

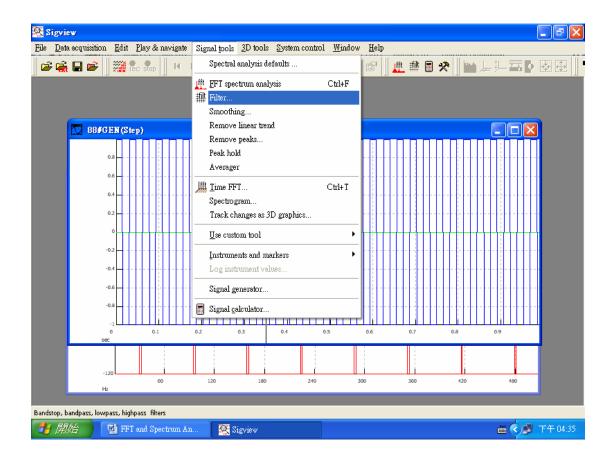

Report:

Discuss the following limitations of the DFT from the results you obtained.

- 1. The signal must be band limited, and the sampling rate must be sufficiently high to avoid aliasing.
- 2. If it necessary to limit the length of the signal for computational purposes, the spectrum will be degraded somewhat by the leakage effect. Leakage is most severe when the simple rectangular window function is used.
- 3. For periodic functions, leakage can be avoided by choosing a proper data length.
- 4. Components lying between discrete frequency lines are subject to error in magnitude due to the "picket-fence" effect.